



InDevR


ENABLING VACCINES & BIOTHERAPEUTICS

From Bottlenecks to Breakthroughs:

Advancing mRNA Vaccine Development
with Same-Day Analytics

Rachel Gao, PhD, Senior Scientist

mRNA Analytical Development Bottlenecks

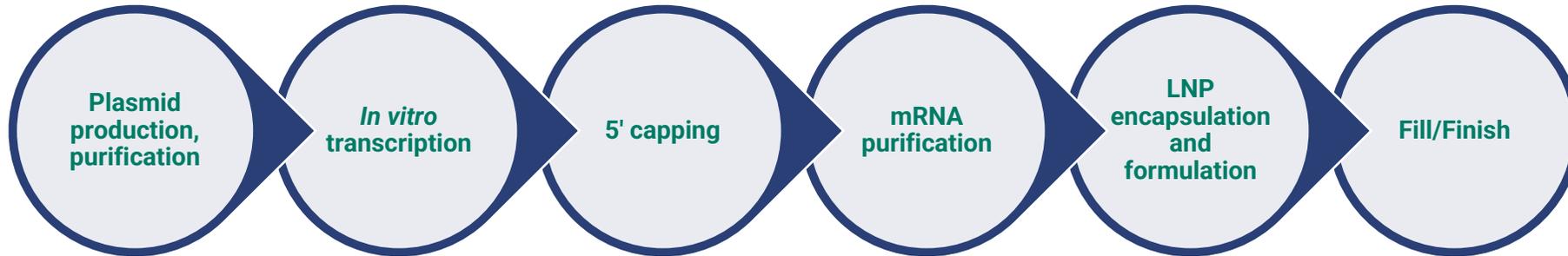

Example Areas of Analytical Technology Constraints

Need

Comprehensive mRNA construct characterization during design, development, and manufacturing

Robust cell-based potency assays

'Validate-able' Release Assays

Challenges

dsRNA, impurities

Method complexity, times to result

Increasingly multivalent formulations

Need Excellent Sensitivity

Increasingly multivalent formulations

Robust, high ease of use, repeatable, GMP-ready

InDevR: Next-Generation Vaccine Analytics



Founded in 2003
Boulder, CO

InDevR's mission is to **eliminate analytical development bottlenecks** in vaccine development and manufacturing

- **2 decades** of streamlining CQA/QC testing for diverse vaccine types (egg-based, cell-based, mRNA)
- **First in industry** to deliver novel strain-responsive off-the-shelf assay kits for **multivalent** influenza antigen testing (HA, NA, NP)
- **Expert Services** division offers tailored solutions for mRNA vaccines as well as traditional vaccines
- Unveiled new cutting-edge facility in Q4 2023 with expanded laboratory capacity

VaxArray Platform Versatility for mRNA Applications



**Multiplexed Assays
Accelerate Analytical
Development Time**



**Standardized
Off-the-Shelf Assay Kits
Empower Reproducibility**



**Custom Assay
Development Services
Available**

Bioprocess

**Rapid multiplexed expressed
protein quantitation**

**Naked mRNA construct identity
and quantity testing**

**Rapid 5'-capping efficiency
assessment**

Drug Product Testing



**Extraction-free LNP-mRNA
identity and quantity testing**

**Rapid multiplexed expressed
protein quantitation**




Clinical/Serology

**Multiplexed quantification of
antibody levels**

Single Instrument, Multiple Assay Options

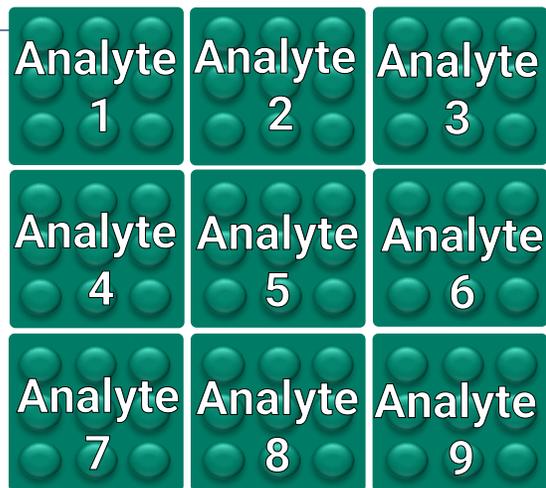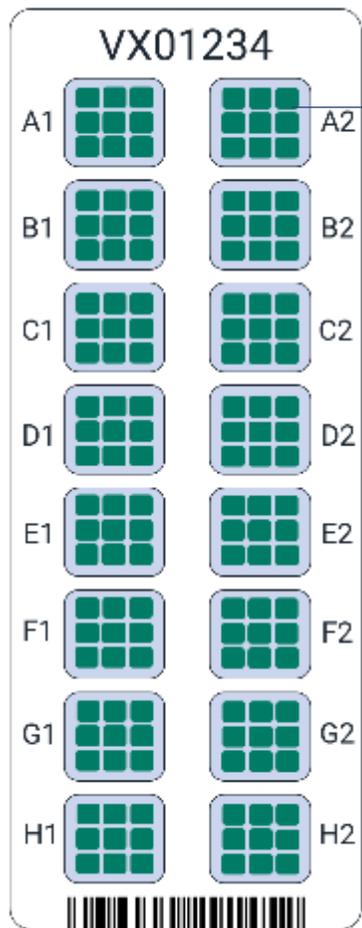

InDevR Expert Services
can tailor custom
assays to your needs,
multiplexing 50+
analytes

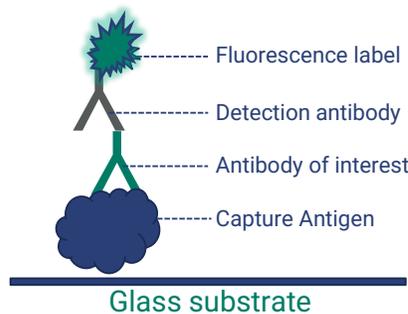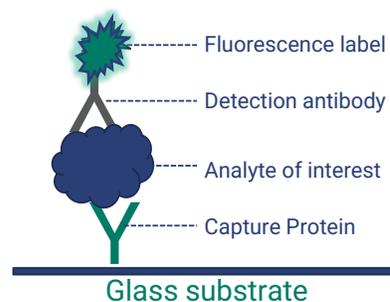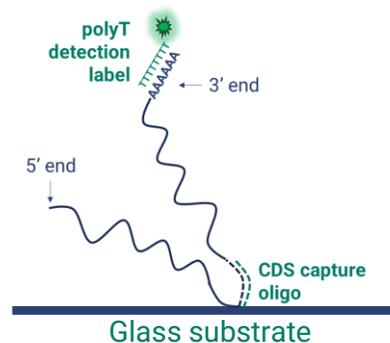

VaxArray mRNA Oligo Assay

- Measure construct-specific mRNA identity and quantity without extraction

VaxArray Antibody Assay

- Measure identity and quantity of protein expression from multiplexed mRNA transfection

VaxArray Antigen Assay



- Measure antibody response from patient serological samples

A close-up photograph of a scientist wearing a white lab coat and blue nitrile gloves. The scientist is holding a pipette with their right hand, positioned over a small vial. The background is a bright, out-of-focus laboratory environment. The overall color palette is dominated by white, blue, and light blue tones.

Applications in mRNA Analytics


In Vitro Potency: Quantify Proteins Post-Transfection

Assess multivalent expressed proteins with high specificity

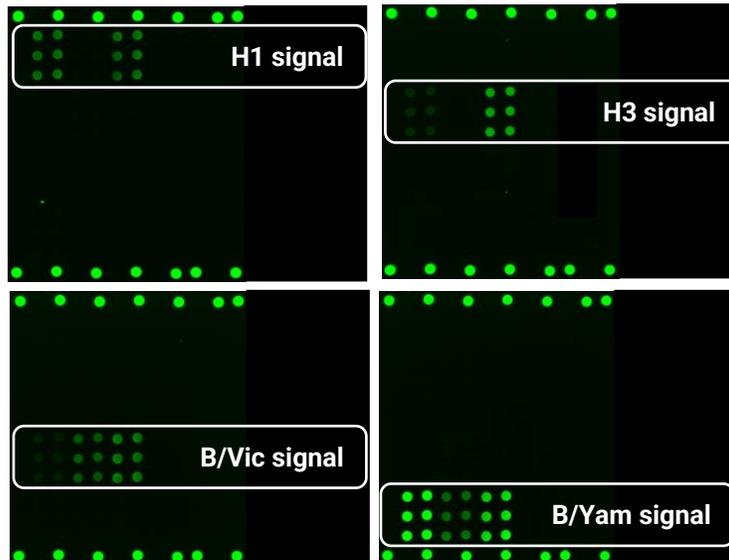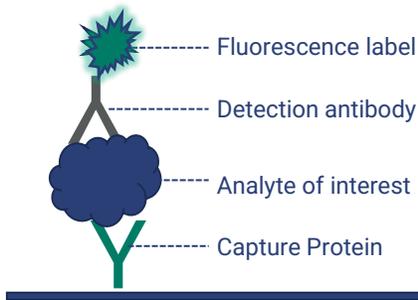


Multiple captures provide reactivity to a broad range of flu strains

Reactivity re-evaluated semi-annually with updated vaccine strains

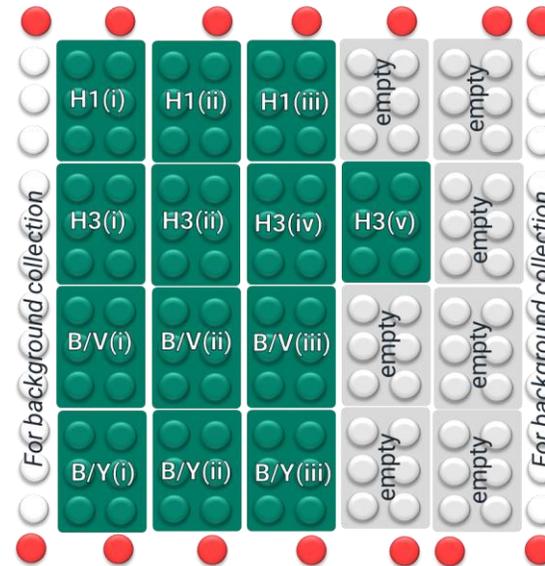

HA microarray layout

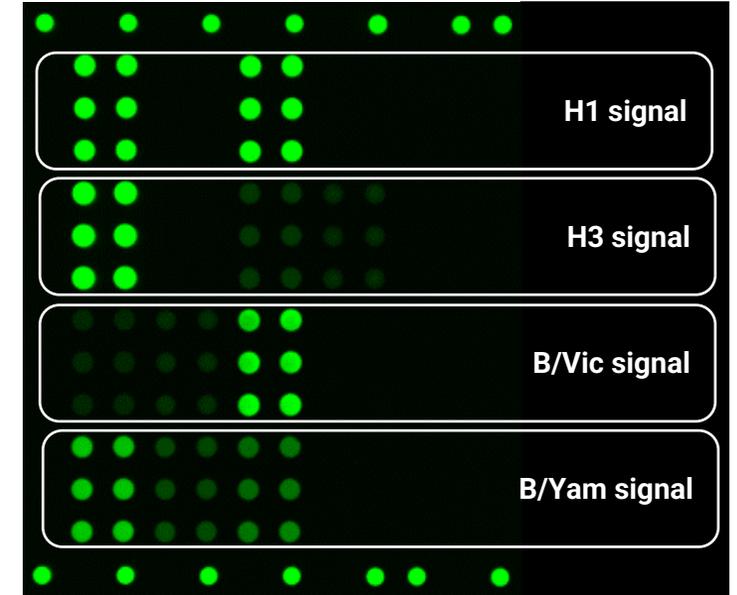


Enables simultaneous quantification of all 4 expressed proteins after quadrivalent mRNA transfection


In Vitro Potency: Quantify Proteins Post-Transfection

Evaluate functional protein during optimization of transfection conditions:

- Input mRNA concentration
 - Sequence optimization
 - Transfection time
 - Batch to batch
-
- HEK293-expressed protein post-transfection with a **4-valent mixture of influenza HA mRNA constructs** measured by VaxArray HA Assay as a function of input mRNA
 - As expected, the different proteins exhibit different expression levels

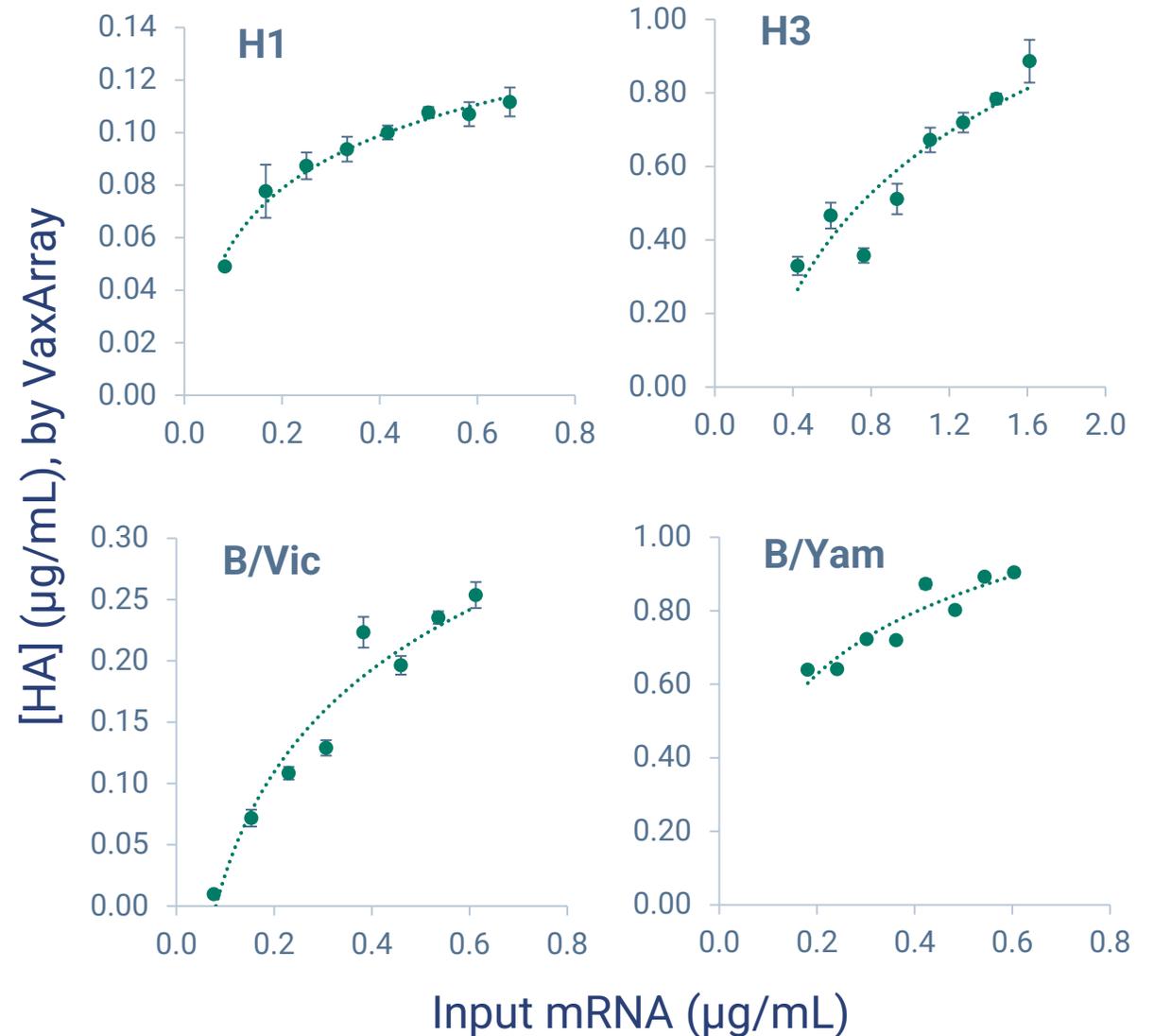

Identity and Quantity of mRNA Constructs

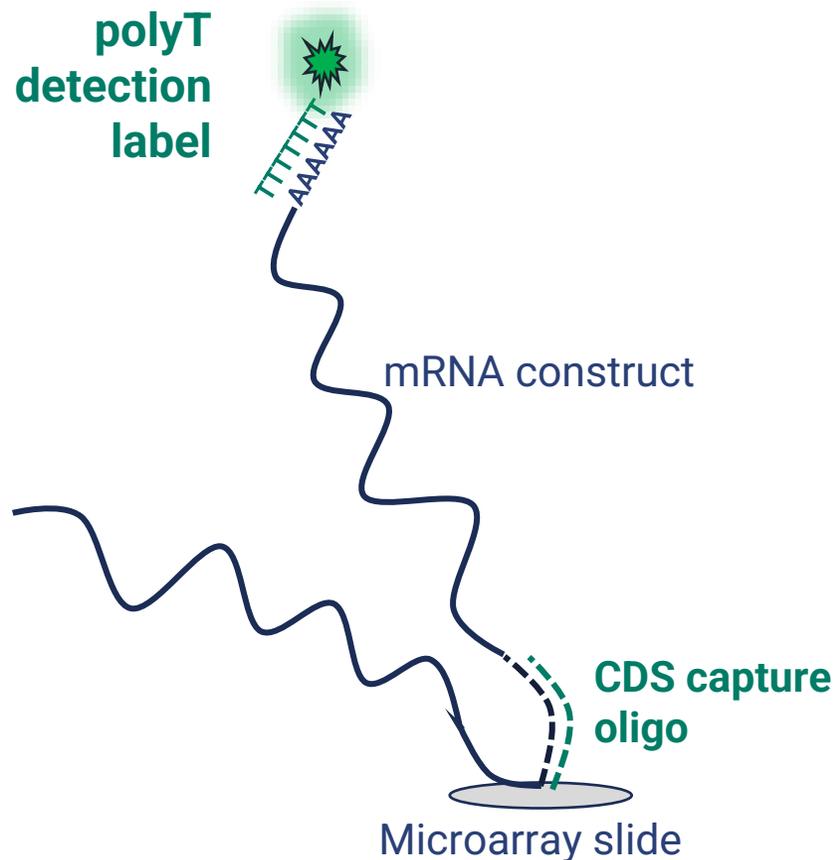

- Identity and quantity of each construct in multivalent mRNA vaccines
- <90-minute time to result
- Capture oligos are designed to target coding regions (CDS) for each construct (ex: each component in a flu vaccine)
- Single detection label for all constructs (custom labels can be designed)
- Works with both naked and LNP-encapsulated mRNA (quick 10-min lysis)
- Can be designed on a custom basis


Identity and Quantity for mRNA Constructs

Case Study: Assay for Identity and quantity in 4-valent flu HA mRNA vaccines


- Oligos designed to capture regardless of strain and codon optimization scheme in a single assay

32 capture oligos, with oligos specific for each vaccine component





- Multiple strains over time spanning **11** vaccine seasons
- 3 unique suppliers (unique UTRs)
- 4 unique codon optimization methods

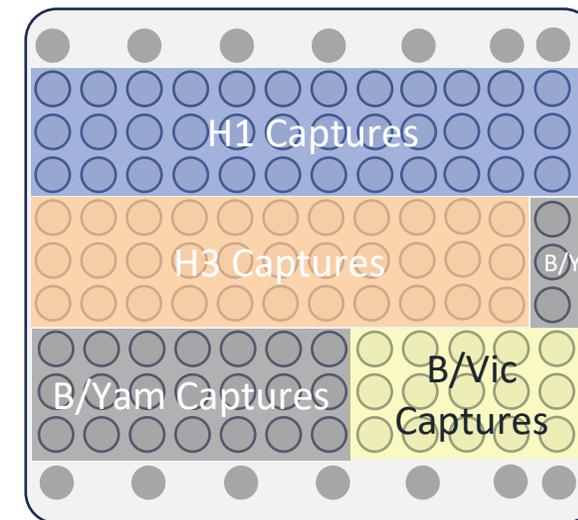

Prototype Array Layout

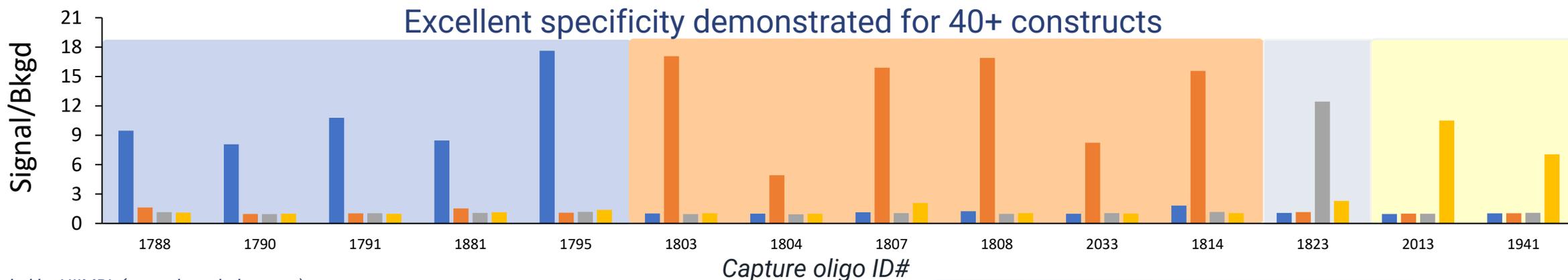


Identity and Quantity for mRNA Constructs

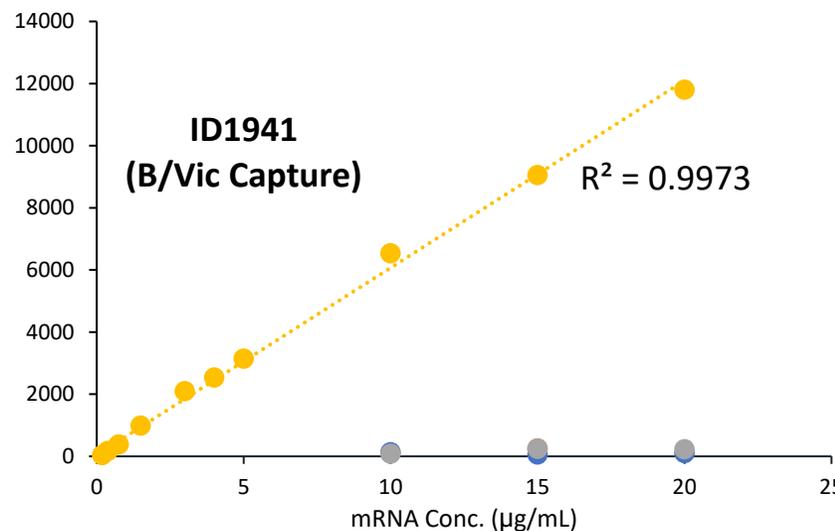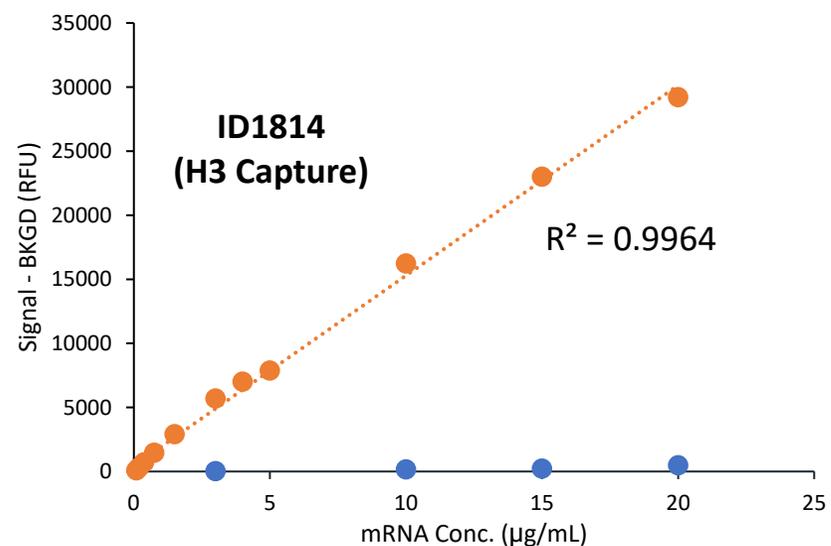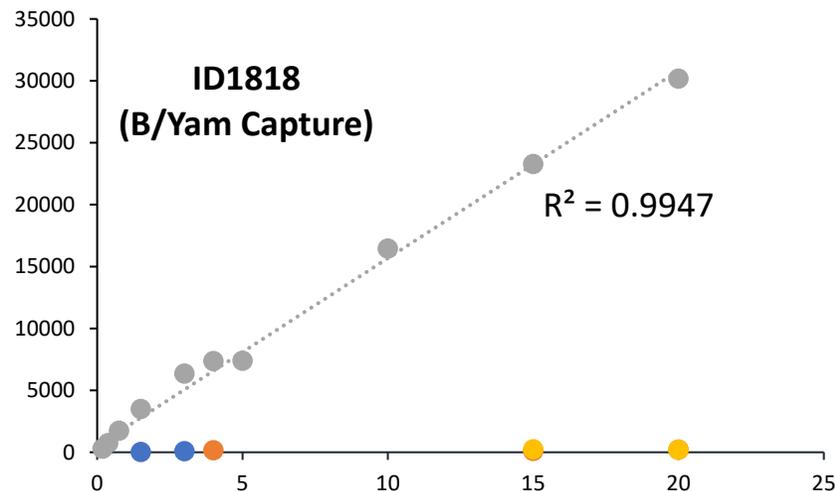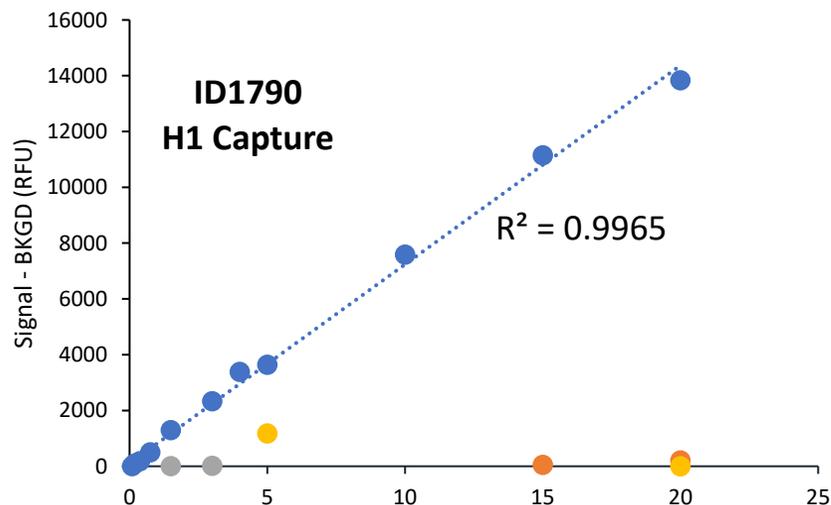

Good linearity over low µg/mL mRNA range

Good specificity: little to no signal on off-target capture oligos


Rapid 5' Capping Assay for Bioprocess Dev't

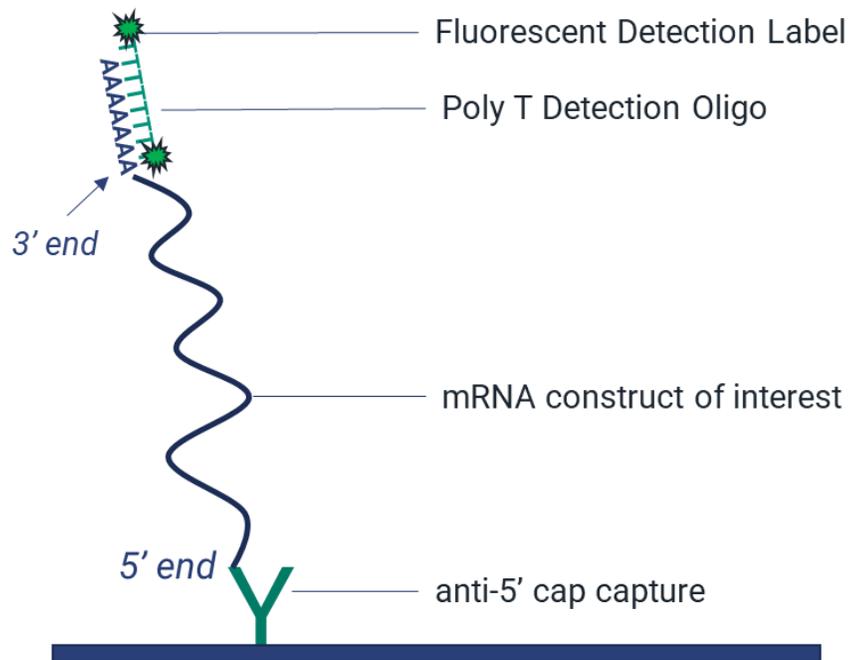

- Capture/detection method is universal to any mRNA construct as long as it has a 5' cap and a 3' polyA tail
- < 2-hour time to result
- Assessment of relative capping or quantitative capping efficiency
- For quantitative analysis, user provides characterized matched standard (same construct) if quantification is desired
- Requires an orthogonal measurement of **total mRNA** on the unknown sample by UV-Vis or other method - therefore material must be reasonably pure

Assay Can Differentiate Wide Range of % Capped

- Contrived samples created by mixing uncapped and capped GFP mRNA construct
- Samples differed by 5% capping; run alongside a dilution series of known capped standard

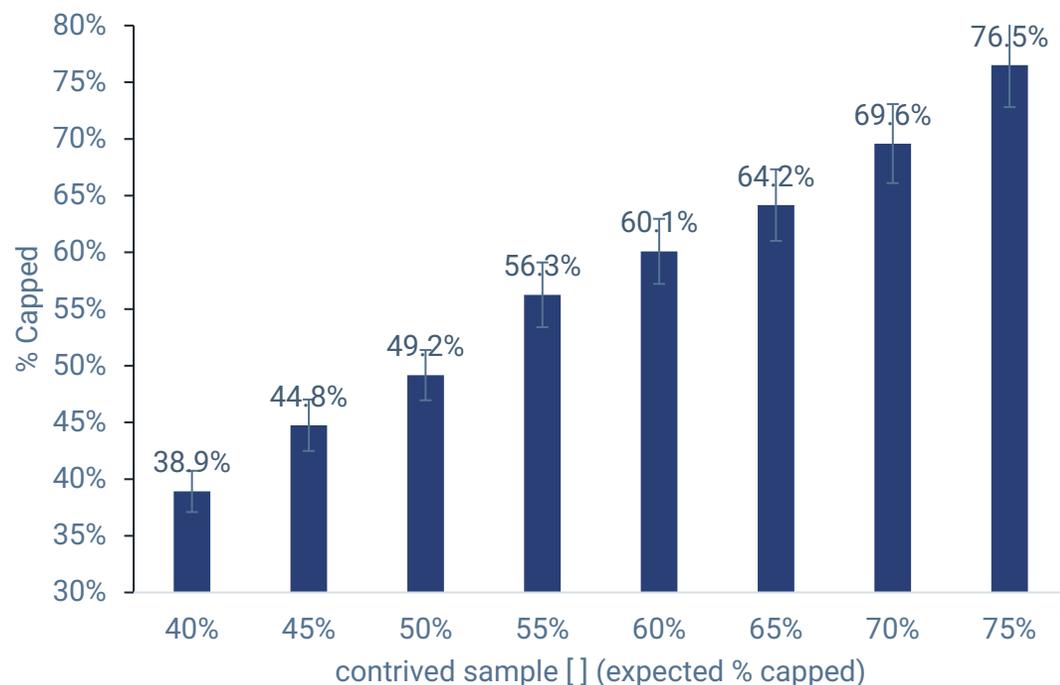

Expected Value	Measured Capping Efficiency (95% CI)	Difference from Expected
40%	38.9 (38.4 – 39.4)%	-1.1%
45%	44.8 (44.1 – 45.3)%	-0.2%
50%	49.2 (48.5 – 49.8)%	-0.8%
55%	56.3 (55.5 – 57.0)%	+1.3%
60%	60.1 (59.3 – 60.8)%	+0.1%
65%	64.2 (63.3 – 65.0)%	-0.8%
70%	69.6 (68.7 - 70.5%)	-0.4%
75%	76.5 (75.5% - 77.5)%	+1.5%

- **All samples statistically differentiated (p-value <0.05 and non-overlapping confidence intervals)**
- **8/8 samples were less than 1.5% difference from expected**



Acknowledgments

Funding (Universal influenza HA mRNA assay): This work was performed under a Project Award Agreement from the National Institute for Innovation in Manufacturing Biopharmaceuticals (NIIMBL) and financial assistance award 70NANB17H002 from the U.S. Department of Commerce, National Institute of Standards and Technology.

Collaborators/Materials Providers

- Dr. Norbert Pardi, University of Pennsylvania, collaborator on NIIMBL funded project
- 5' capping assay, materials provided and used during development:
 - North Carolina State University/BTEC, Dr. Jennifer Pancorbo
 - Cisterna, Dr. Hari Bhaskaran
 - Aldevron

InDevR's Expert Services and R&D Teams



We Want to Hear From You



Scan to connect with InDevR's scientists

